



Ultra Low Temperature RFID Tag Application Guide





Ultra Low Temperature RFID Tag Application Guide

Supplies Required:

- 9x12 Poly Bag
- 12x12 Poly Bag
- Tag Flag
- RFID Tag
- Stapler (Optional)



Process:

Step 1. Locate product and determine bag size needed based on product size. (9"x12" or 12"x12")


Step 3. Associate product and place inside bag.

Step 4. Take tag flag and a RFID tag. On the tag flag locate the paper covering at the top of the tag flag.





Ultra Low Temperature RFID Tag Application Guide

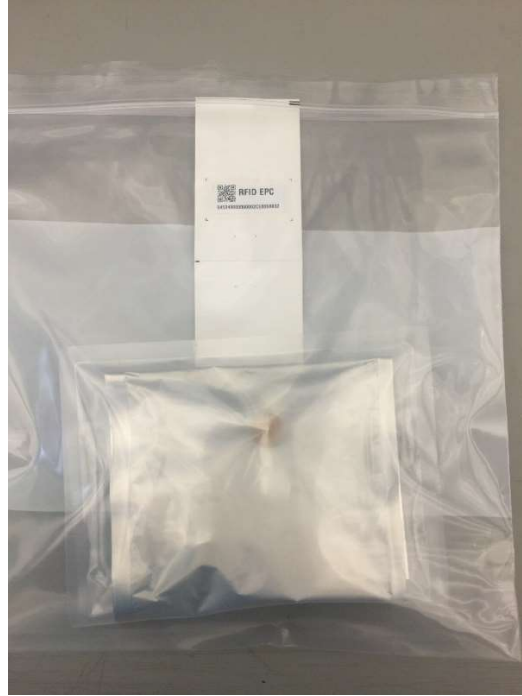

Step 7. Place inside freezer blue bin with the RFID tag and product label facing outward. (Optional: Staple the flag to the bag to keep it upright. Avoid stapling the RFID tag itself)

Step 8. Ensure products are facing outward and that the plastic flags are staggered and not stacked one in front of another.